



HOW TO CREATE A NEW USER IN PARAVISION 6

1. Login as nmrsu
2. In the menu select *Applications* → *Administration* → *User and Groups*

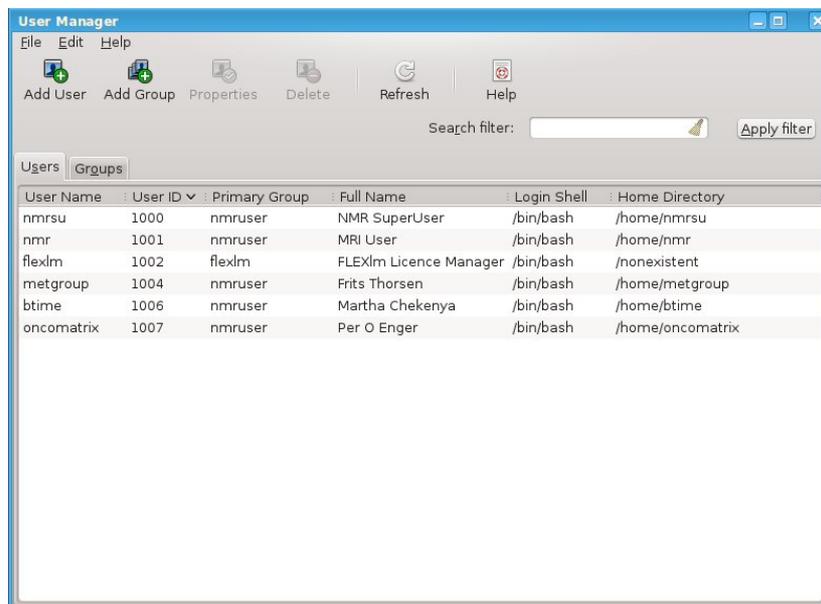


3. Press *Add User*

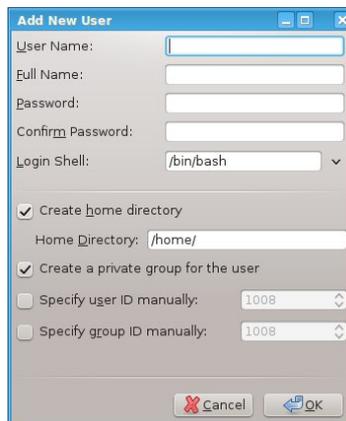

4. Select a *User Name*
 - Since we are creating one user name per research group, the User Name should preferably be the name of the group.

5. Write the *Full Name* of the groups' PI.
6. Choose a password that is easy for the group to remember
7. Press OK
8. Click on *Properties*
9. Make the user be part of the *nmruser* group by clicking on *nmruser*
10. Make the user be part of the *nmrsu* group by clicking on *nmrsu*
11. Make *nmruser* be the primary group of the user (this is required for PV to work).
12. Close the user *Manager*
13. Logout nmrsu
14. Login as the new user
15. To start PV for the first time:
 - Open a shell/terminal
 - `cd /opt/PV6.01./pv`
 - Select Y for all the questions
 - ParaVision is now starting

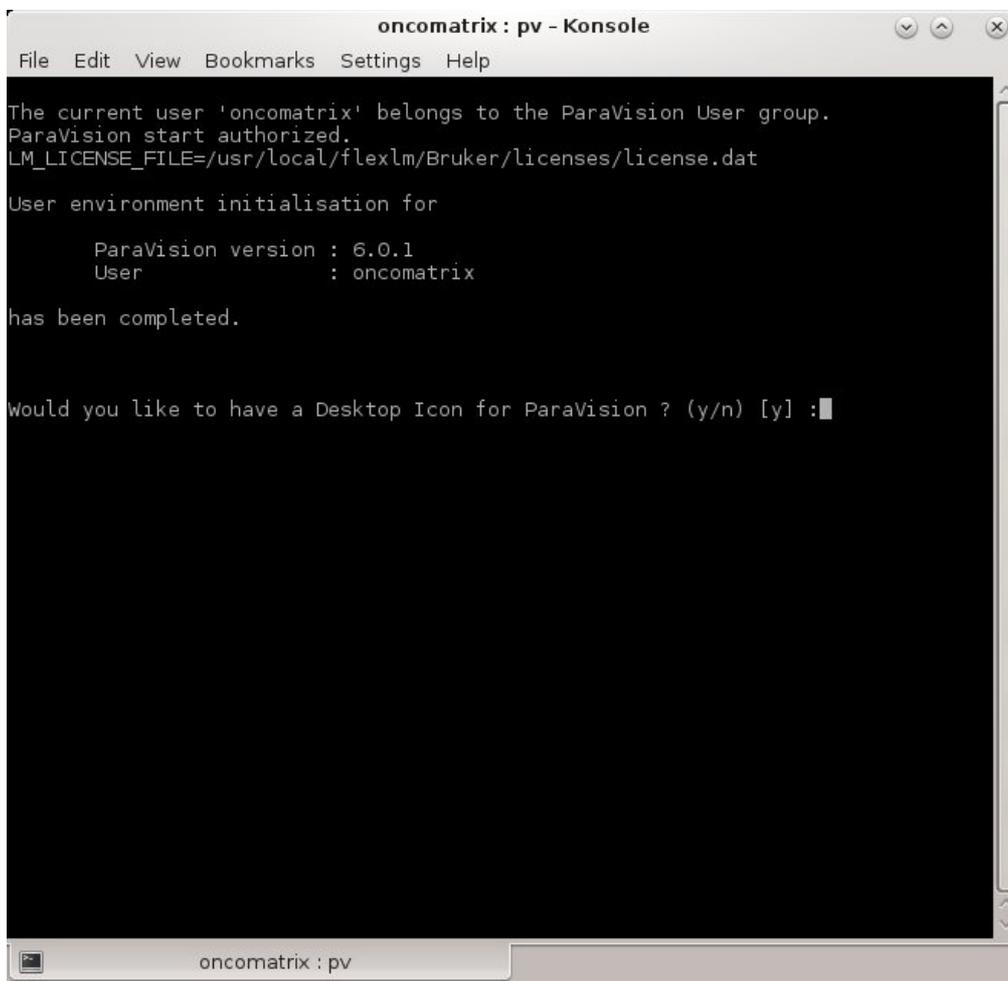

```
oncomatrix : pv - Konsole
File Edit View Bookmarks Settings Help

The current user 'oncomatrix' belongs to the ParaVision User group.
ParaVision start authorized.
LM_LICENSE_FILE=/usr/local/flexlm/Bruker/licenses/license.dat

User environment initialisation for

    ParaVision version : 6.0.1
    User                : oncomatrix

has been completed.

Would you like to have a Desktop Icon for ParaVision ? (y/n) [y] :
```

16. Close ParaVision
17. Start ParaVision again by clicking on the PV icon on the Desktop
18. Run a Tripilot to see that everything is ok.