



## CHECKLIST FOR WHEN YOU ARE DONE SCANNING

<p><b>Log into the blue book in the control room</b> Record date and time of scanning, project title, level of liquid N<sub>2</sub> and He, user name</p>
<p><b>Turn off power in the first cabinet with high power electronics</b> FIRST, press the top (on/off) button, under local (h.v) THEN, switch the system off using the switch above power</p>
<p><b>Turn off the hot water circulation system</b> Make sure the heating blanket remains connected to the tubing</p>
<p><b>Disconnect all anesthesia gases and Turn off isoflurane on the anesthesia vaporizer</b> Applies to the MR and animal preparation room</p>
<p><b>Disconnect the battery from the temperature module and connect it to the charging port on the respiration module</b></p>
<p><b>Clean the animal bed and table</b> Spray with liquid in the red bottle, leave acting for 10 min, rinse with water in the blue bottle</p>
<p><b>Remove and store any equipment that is particular to your experiment</b> Use cabinets in the storage room next to the animal preparation room</p>
<p><b>Tidy up behind you</b> Remove all syringes, catheters, contrast agents, etc...</p>
<p><b>Switch off the light in the MR room</b> Light source in the MR room heats up the room</p>
<p><b>Lock the doors behind you</b> Lock the main door the MR facility</p>